

UNDECAGOLD



95 Horse Block Road, Yaphank NY 11980-9710
Tel: (877) 447-6266 (Toll-Free in US) or (631) 205-9490 Fax: (631) 205-9493
Tech Support: (631) 205-9492 tech@nanoprobes.com
www.nanoprobes.com

PRODUCT INFORMATION

MONOAMINO UNDECAGOLD LABELING REAGENT

Product Name: MONOAMINO UNDECAGOLD
Catalog Number: 2031
Appearance: Orange-brown solid
Revision: 1.3 (March 2000)

Congratulations on your acquisition of a revolutionary new gold immunoreagent: the MONOAMINO UNDECAGOLD labeling kit. With this reagent you can label glycoproteins, or any other biomolecule containing a suitable reactive group, with UNDECAGOLD. Because UNDECAGOLD is a discrete molecular compound and not a colloidal gold preparation, conjugates prepared with this reagent have several advantages over colloidal gold conjugates (see below).

CONTENTS

Product Information
Product Specifications
Thiol Caution
Labeling Glycoproteins with UNDECAGOLD
Labeling Other Molecules with UNDECAGOLD
General Considerations for Staining with UNDECAGOLD Reagents
Using Stains with UNDECAGOLD
Electron Microscopy Labeling with UNDECAGOLD
 1. Cells in Suspension
 2. Thin Sections
Special Considerations for Viewing UNDECAGOLD in the Electron Microscope
Silver Enhancement of UNDECAGOLD for EM
Light Microscopy with UNDECAGOLD
Immunoblotting with UNDECAGOLD
References

Warning: For research use only. Not recommended or intended for diagnosis of disease in humans or animals. Do not use internally or externally in humans or animals. Non radioactive and non carcinogenic.

PRODUCT INFORMATION

UNDECAGOLD is the smallest gold label available, prepared using a discrete gold compound rather than a colloid.¹ This kit contains the UNDECAGOLD particle with a single primary amine functionality incorporated into a ligand on the surface of the gold particle; this has a specific reactivity towards aldehyde groups, with which it forms Schiff bases. These may then be reduced to amines. This

process may be used to label carbohydrate moieties in glycoproteins, as shown on the next page (figure 1). It may also be used with cross-linking reagents to label other groups, such as primary amines. The reagent as supplied has been lyophilized from methanol; dissolution in the reaction solvent or buffer will produce a solution of activated UNDECAGOLD. UNDECAGOLD conjugates may be used in electron microscopy and some immunoblotting and light microscopy applications. They are stable to wide ranges of pH and ionic strength, and are not radioactive or carcinogenic. This product should be stored at -20°C.

PRODUCT SPECIFICATIONS

MONOAMINO UNDECAGOLD is supplied lyophilized from methanol solution. The dark brown solid may be dissolved in aqueous buffer systems such as phosphate buffered saline or ammonium acetate, and is also soluble in alcohols, acetone, DMSO, dichloromethane and other organic solvents. Extinction coefficients at specific wavelengths are given below for methanol solution:

<u>WAVELENGTH (nm)</u>	<u>EXTINCTION COEFFICIENT*</u>
280	1.677 X 10 ⁵
420	0.471 X 10 ⁵

*Measured for 5 X 10⁻⁶ M solution in methanol.

50 nmol of the product is supplied. The amounts required will vary according to the system under study, but in general a 5-fold to 10-fold excess of UNDECAGOLD is recommended; the reagent supplied will therefore label up to 0.5 - 1.0 mg of a molecule with one labeling site and a molecular weight of 100 000.

THIOL CAUTION

UNDECAGOLD particles degrade upon exposure to concentrated thiols such as β-mercaptoethanol or dithiothreitol. If such reagents must be used, concentrations should be kept below 1 mM and exposure restricted to 10 minutes or less.

LABELING GLYCOPROTEINS WITH UNDECAGOLD

MONOAMINE UNDECAGOLD is reactive towards aldehydes. It may be used to label carbohydrate moieties of glycoproteins,² as shown in figure 1 (next page).

Sufficient UNDECAGOLD reagent is supplied to label up to 10 nmol of aldehyde groups (for example, 1.0 mg of a compound with one aldehyde and a molecular weight of 100,000). UNDECAGOLD is stable for several weeks in most aqueous buffer solutions. For purification and isolation steps alternative buffers may be substituted for those given; however, the labeling reaction itself should be performed using the buffers and conditions specified.

To assist structural preservation, it is usually helpful to fix proteins before labeling. If glutaraldehyde or other aldehyde-containing fixing reagents are used, these should be quenched before labeling. This may be achieved by reacting with sodium borohydride (15-20 mole equivalents per glutaraldehyde) at 4°C for 1.5 h.² Alternatively, incubate the specimens for 5 minutes in 50 mM glycine solution in PBS (pH 7.4), or treat with ammonium chloride (50 mM).

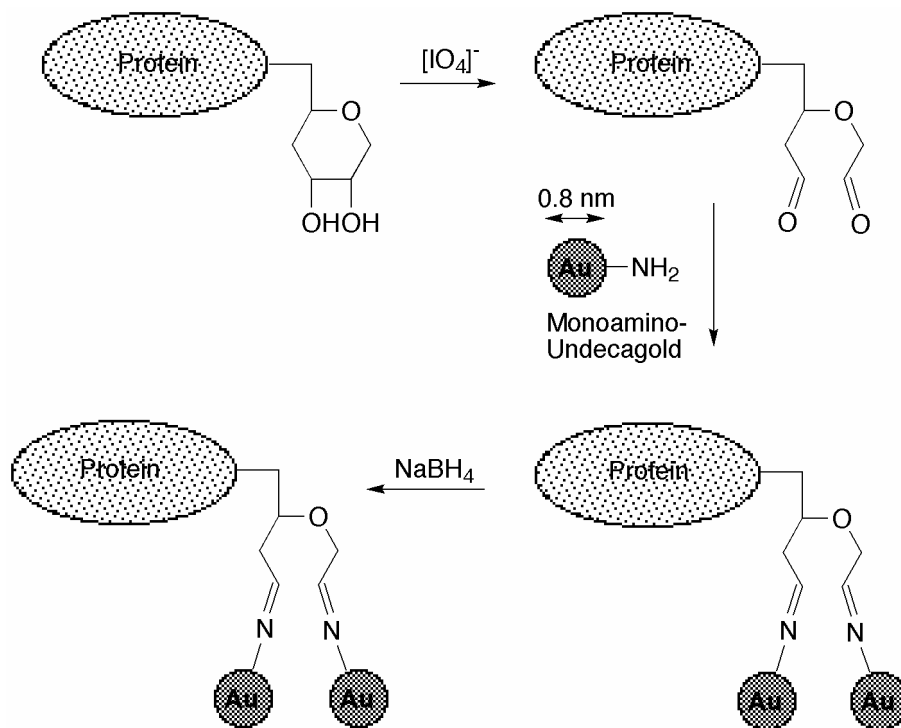


Figure 1: Labeling of a carbohydrate moiety with MONOAMINO UNDECAGOLD.²

1. If the labeling site is in the form of a carbohydrate group, it must be oxidized to produce aldehydes. Incubate specimen for 1.5 h at 4°C with 1000-2000 mole equivalents of sodium periodate per carbohydrate in 20 mM PIPES.Na.
2. Isolate oxidized glycoprotein by dialysis or gel filtration chromatography to remove sodium periodate. Use a gel such as Amicon GH-25, which has an exclusion cut-off at molecular weight 3,000. As the eluent, use 20mM PIPES.Na, with 150 mM sodium chloride. The oxidized glycoprotein will be eluted in the void volume as the first sharp peak in the trace. Combine the fractions containing protein; the total amount may be calculated from the optical density.
3. Dissolve the UNDECAGOLD reagent in 1 ml buffer. Sufficient reagent is supplied to label up to 10 nmol of aldehydes; if you are using a smaller amount, use a proportionately smaller amount of the NANOGOLD™ reagent.
4. Add the MONOAMINO UNDECAGOLD solution to the oxidized glycoprotein (6 to 10 mole equivalents per carbohydrate chain). Incubate for 1 hour at 4°C.
5. Reduce the Schiff base linkages with excess NaBH₄; use 500-1000 mole equivalents per carbohydrate chain, and allow to react for 30 minutes at 4°C; then quench with excess acetone.
6. Separate the unbound gold particles from the labeled glycoprotein using gel exclusion chromatography. The UNDECAGOLD conjugate may be effectively isolated using a medium such as Pharmacia Superose 6 or 12 (which fractionate a wide range of molecular weights) or Amicon GCL-90 (which excludes molecules of mass 30,000 or greater). Concentrate the reaction mixture to a suitably small volume using membrane centrifugation (e.g. Amicon Centricon-30 system). Elute with 0.25 M ammonium acetate (nominal pH 7.7); monitor at 280 nm. The first, faintly colored peak is the conjugate, while the second, darker band is unbound UNDECAGOLD particles. For even higher purity, repeat this process one time.

UNDECAGOLD conjugates should be stored in 0.02 M phosphate buffer with 150 mM sodium chloride, or other buffer solutions usually used with the protein under study. If they are to be stored longer than three days, add 0.1 % bovine serum albumin and 0.05 % sodium azide to prevent bacterial contamination and to prevent the protein from adhering to the surfaces of the storage vessel.

If the molecule to be labeled already contains aldehydes, the procedure above should be followed without the periodate oxidation and removal.

LABELING OTHER MOLECULES WITH UNDECAGOLD

MONOAMINO UNDECAGOLD may be used to label other groups if it is used with a suitable cross-linker. As an example, the use of bis (sulfosuccinimidyl) suberate (BS³) as a cross linker to label a protein with a primary amine is described below.³ The reaction works best in concentrated protein solutions (i.e. smaller reaction volumes).

1. Dissolve the protein in 0.02 M sodium phosphate with 150 mM sodium chloride at pH 7.4 (0.5 ml).
2. Dissolve the UNDECAGOLD reagent in 1 ml buffer. Sufficient reagent is supplied to label up to 10 nmol of sites; if you are using a smaller amount, use a proportionately smaller amount of the UNDECAGOLD reagent.
3. Add the MONOAMINO UNDECAGOLD solution to the protein (6 to 10 mole equivalents per mole equivalent amine labeling site). Mix thoroughly.
4. Dissolve BS³ in DMSO (0.1 ml) and dilute to 2 ml with 0.02 M phosphate buffer with 150 mM NaCl, pH 7.4. Use sufficient BS³ to give a concentration of 1-2 mM in the reaction mixture (0.5 mg). Once dissolved, the cross-linking reagent should be used immediately. Hydrolysis will occur within a few hours.
5. Add the BS solution to the protein/UNDECAGOLD solution, and incubate for 10 mins at room temperature.
6. Quench with excess 10 mM tris buffer at pH 7.0.
7. Separate the unbound gold particles from the labeled glycoprotein using gel exclusion chromatography. The UNDECAGOLD conjugate may be effectively isolated using a medium such as Pharmacia Superose 6 or 12 (which fractionate a wide range of molecular weights) or Amicon GCL-90 (which excludes molecules of mass 30,000 or greater). Concentrate the reaction mixture to a suitably small volume using membrane centrifugation (e.g. Amicon Centricon-30 system). Elute with 0.25 M ammonium acetate (nominal pH 7.7); monitor at 280 nm. The first, faintly colored peak is the conjugate, while the second, darker band is unbound UNDECAGOLD particles. For even higher purity, repeat this process one time.

Other cross-linking reagents may be used to link the MONOAMINO UNDECAGOLD to other functionalities. The buffer and reagent concentrations used for this reaction, and the incubation time, may be varied according to the nature of the molecule under study.

GENERAL CONSIDERATIONS FOR IMMUNOSTAINING WITH UNDECAGOLD REAGENTS

Basically, normal methodologies may be used successfully with UNDECAGOLD immunoreagents. The concentration of antibody and gold is similar to other commercial preparations of colloidal gold antibodies. Therefore similar dilutions and blocking agents are appropriate.

The major difference will be in the results:

UNDECAGOLD is an extremely uniform 0.8 nm diameter gold particle ($\pm 10\%$).

UNDECAGOLD conjugates contain absolutely no aggregates. This is in sharp contrast to other colloidal gold conjugates that usually are prepared by centrifugation to remove the largest aggregates and frequently contain smaller aggregates.

Close to 1 UNDECAGOLD particle to 1 Fab (or IgG) make this product distinct from the 0.2 - 10 variable stoichiometry of other colloidal gold antibody preparations.

UNDECAGOLD particles do not have affinity to proteins as do other other colloidal golds. This reduces background and false labeling.

USING STAINS WITH UNDECAGOLD

Because the 0.8 nm UNDECAGOLD particles are so small, over staining with OsO₄, uranyl acetate or lead citrate will obscure direct visualization of individual UNDECAGOLD particles, and therefore these stains should not be used. Only light staining with a low atomic number stain, such as NANOVANTM, a Vanadium based negative stain, should be used.

ELECTRON MICROSCOPY IMMUNOLABELING WITH UNDECAGOLD

If aldehyde-containing reagents have been used for fixation, these must be quenched before labeling. This may be achieved by incubating the specimens for 5 minutes in 50 mM glycine solution in PBS (pH 7.4). Ammonium chloride (50 mM) or sodium borohydride (0.5 - 1 mg/ml) in PBS may be used instead of glycine.

Cells in Suspension

1. Optional fixing of cells: e.g., with glutaraldehyde (0.05 - 1% for 15 minutes) in PBS. Do not use Tris buffer since this contains an amine. After fixation, centrifuge cells (e.g. 1 ml at 10^7 cells/ml) at 300 X g, 5 minutes; discard supernatant; resuspend in 1 ml buffer. Repeat this washing (centrifugation and resuspension) 2 times.
2. Incubate cells with 0.02 M glycine in PBS (5 mins). Centrifuge, then resuspend cells in PBS-BSA buffer (specified below) for 5 minutes.
4. Wash cells using PBS-BSA as described in step 1 (2 X 5 mins). Resuspend in 1 ml Buffer 1.
3. Place 50 - 200 μ l of cells into Eppendorf tube. Dilute UNDECAGOLD conjugate ~ 50 times in PBS-BSA buffer and add 30 μ l to cells; incubate for 30 minutes with occasional shaking (do not create bubbles which will denature proteins).
5. Wash cells in PBS-BSA as described in step 1 (2 X 5 mins).
6. Fix cells and antibodies using a final concentration of 1% glutaraldehyde in PBS for 15 minutes. Then remove fixative by washing with buffer 1 (3 X 5 mins).

PBS-BSA Buffer:

20 mM phosphate
150 mM NaCl
pH 7.4
0.5% BSA
0.1% gelatin (high purity)

Optional, may reduce background:

0.5 M NaCl
0.05% Tween 20

PBS Buffer:

20 mM phosphate
150 mM NaCl
pH 7.4

Negative staining may be used for electron microscopy of small structures or single molecules which are not embedded. Negative stain must be applied after the silver enhancement. NANOVAN™ negative stain is specially formulated for use with UNDECAGOLD reagents; it is based on vanadium, which gives a lighter stain than uranium, lead or tungsten-based negative stains and allows easier visualization of UNDECAGOLD particles with little or no silver enhancement.

Thin Sections

Labeling with UNDECAGOLD may be performed before or after embedding.⁴ Labeling before embedding and sectioning (the pre-embedding method) is used for the study of surface antigens, particularly small organisms such as viruses budding from host cells. It gives good preservation of cellular structure, and subsequent staining usually produces high contrast for study of the cellular details. Labeling after embedding and sectioning (the post-embedding method) allows the antibody access to the interior of the cells, and is used to label both exterior and interior features. The procedures for both methods are described below.

Thin sections mounted on grids are floated on drops of solutions on parafilm or in well plates. Hydrophobic resins usually require pre-etching.

PROCEDURE FOR PRE-EMBEDDING METHOD:⁴

1. Float on a drop of water for 5 - 10 minutes.
2. Incubate cells with 1 % bovine serum albumin in PBS buffer at pH 7.4 for 5 minutes; this blocks any non-specific protein binding sites and minimizes non-specific antibody binding.
3. Rinse with PBS-BSA (1 min).

4. Incubate with UNDECAGOLD conjugate diluted 1/40 - 1/200 in PBS-BSA with 1 % normal serum from the same species as the UNDECAGOLD labeled antibody, for 10 minutes to 1 hour at room temperature.
5. Rinse with PBS-BSA (3 X 1 min), then PBS (3 X 1 min).
6. Postfix with 1 % glutaraldehyde in PBS (10 mins).
7. Rinse in deionized water (2 X 5 min).
8. Dehydrate and embed according to usual procedure. Use of a low-temperature resin (eg. Lowicryl) is recommended.
9. Stain (uranyl acetate, lead citrate or other positive staining reagent) as usual before examination.

Silver enhancement may be performed before or after embedding (see below); it should be completed before postfixing or staining with osmium tetroxide, uranyl acetate or similar reagents is carried out.

PROCEDURE FOR POST-EMBEDDING METHOD:⁴

1. Prepare sections on plastic or carbon-coated nickel grid. Float on a drop of water for 5 - 10 minutes.
2. Incubate with 1 % solution of bovine serum albumin in PBS buffer at pH 7.4 for 5 minutes to block non-specific protein binding sites.
3. Rinse with PBS-BSA (1 min).
4. Incubate with UNDECAGOLD conjugate diluted 1/40 - 1/200 in PBS-BSA with 1 % normal serum from the same species as the UNDECAGOLD labeled antibody, for 10 minutes to 1 hour at room temperature.
5. Rinse with PBS (3 X 1 min).
6. Postfix with 1 % glutaraldehyde in PBS at room temperature (3 mins).
7. Rinse in deionized water for (2 X 5 min).
8. If desired, contrast sections with uranyl acetate and/or lead citrate before examination.

Silver enhancement may also be used to render the UNDECAGOLD particles more easily visible (see below), especially if stains such as uranyl acetate or lead citrate are applied. If used, it should be completed before these stains are applied.

PBS-BSA Buffer:

20 mM phosphate
150 mM NaCl
pH 7.4
0.5% BSA
0.1% gelatin (high purity)

Optional, may reduce background:

0.5 M NaCl
0.05% Tween 20

PBS Buffer:

20 mM phosphate
150 mM NaCl
pH 7.4

SPECIAL CONSIDERATIONS FOR VIEWING UNDECAGOLD IN THE ELECTRON MICROSCOPE

UNDECAGOLD is the smallest gold probe commercially available, being just 0.8 nm in diameter. A high resolution instrument such as a Scanning Transmission Electron Microscope (STEM) is required for visualization; in a conventional TEM the UNDECAGOLD particles are not visible. With careful work, however, UNDECAGOLD may be seen directly in the STEM. However, achieving the high resolution necessary for this work may require new demands on your equipment and technique. Several suggestions follow:

1. Before you start a project with UNDECAGOLD it is helpful to see it so you know what to look for. Dilute the UNDECAGOLD stock 1:5 in methanol and apply 4 μ l to a grid for 1 minute. Allow to dry.
2. View UNDECAGOLD using a full width scan of 128 nm or less; this will give sufficient magnification for visualization.
3. UNDECAGOLD is sensitive to beam damage (contrary to NANOGOLD[®] which is very beam-resistant); the behavior of UNDECAGOLD in the STEM has been described in the literature.² Image at approximately 200 e⁻ \AA^{-2} .

4. In order to operate at high magnification, thin carbon film over fenestrated holey film is recommended. Many plastic supports are unstable under these conditions of high magnification/high beam current and carbon is therefore preferred. Contrast is best using thinner films.

SILVER ENHANCEMENT OF UNDECAGOLD FOR EM

UNDECAGOLD will nucleate silver deposition resulting in a dense particle 2-20 nm in size or larger depending on development time. However, silver enhancement will be slower and much less uniform than with larger gold particles such as NANOGOLD®. If specimens are to be embedded, silver enhancement is usually performed after embedding, although it may be done first. It must be completed before any staining reagents such as osmium tetroxide, lead citrate or uranyl acetate are applied, since these will nucleate silver deposition in the same manner as gold and produce non-specific staining.

Our LI SILVER silver enhancement system is convenient and not light sensitive, and suitable for all applications. Improved results in the EM may be obtained using HQ SILVER, which is formulated to give slower, more controllable particle growth and uniform particle size distribution.

Specimens must be thoroughly rinsed with deionized water before silver enhancement reagents are applied. This is because the buffers used for antibody incubations and washes contain chloride ions and other anions which form insoluble precipitates with silver. These are often light-sensitive and will give non-specific staining. To prepare the developer, mix equal amounts of the enhancer and initiator immediately before use. UNDECAGOLD will nucleate silver deposition resulting in a dense particle 2-20 nm in size or larger depending on development time. Use nickel grids (not copper).

The procedure for immunolabeling should be followed up to step 6 as described above. Silver enhancement is then performed as follows:

1. Rinse with deionized water (2 X 5 mins).
2. Float grid with specimen on freshly mixed developer for 1-4 minutes, or as directed in the instructions for the silver reagent. More or less time can be used to control particle size. A series of different development times should be tried, to find the optimum time for your experiment.
3. Rinse with deionized water (3 X 1 min).
4. Mount as usual.

LIGHT MICROSCOPY WITH UNDECAGOLD

For most light microscopy applications we recommend NANOGOLD® conjugates, since these usually give more intense, specific signals. UNDECAGOLD must be developed with a silver enhancement reagent before it is visible in the light microscope. Our LI SILVER is convenient for this as it is not light sensitive and the degree of enhancement may be monitored readily.

Samples must be rinsed with deionized water before silver enhancement. This is because the reagent contains silver ions in solution, which react to form a precipitate with chloride, phosphate and other anions which are components of buffer solutions. The procedure for immunolabeling with UNDECAGOLD and silver enhancement is given below.

1. Spin cells onto slides using Cytospin, or use paraffin section.
2. Incubate with 1 % solution of bovine serum albumin in PBS (PBS-BSA) for 10 minutes to block non-specific protein binding sites.
3. Incubate with primary antibody, diluted at usual working concentration in PBS-BSA (1 hour or usual time)
4. Rinse with PBS-BSA (3 X 2 min).
5. Incubate with UNDECAGOLD reagent diluted 1/40 - 1/200 in PBS-BSA with 1 % normal serum from the same species as the UNDECAGOLD reagent, for 1 hour at room temperature.
6. Rinse with PBS (3 X 5 min).
7. Postfix with 1 % glutaraldehyde in PBS at room temperature (3 mins).
8. Rinse with deionized water (3 X 1 min).
9. Develop specimen with freshly mixed developer for 5-20 minutes, or as directed in the instructions for the silver reagent. More or less time can be used to control intensity of signal. A series of different development times may be used, to find

the optimum enhancement for your experiment; generally a shorter antibody incubation time will require a longer silver development time.

10. Rinse with deionized water (2 X 5 mins).
11. The specimen may now be stained if desired before examination, with usual reagents.

PBS-BSA Buffer:

20 mM phosphate
150 mM NaCl
pH 7.4
0.5% BSA
0.1% gelatin (high purity)

Optional, may reduce background:

0.5 M NaCl
0.05% Tween 20

PBS Buffer:

20 mM phosphate
150 mM NaCl
pH 7.4

To obtain an especially dark silver signal, the silver enhancement may be repeated with a freshly mixed portion of developer.

IMMUNOBLOTTING

For immunoblotting we recommend our NANOGOLD[®] conjugates, since these give more sensitive detection than UNDECAGOLD conjugates with a higher signal-to-noise ratio upon silver enhancement. The basic procedure for gold immunoblotting has been described by Moeremans et al⁵, which may be followed. For best results, the membrane should be hydrated before use by simmering in gently boiling water for 15 minutes. Best results are obtained when the antigen is applied using a 1 µl capillary tube. A suitable procedure for immunoblots is as follows:

1. Spot 1 µl dilutions of the antigen in buffer 4 onto hydrated nitrocellulose membrane. Use an antigen concentration range from 100 to 0.01 pg / µl.
2. Block with buffer 1 for 30 minutes at 37°C.
3. Incubate with primary antibody according to usual procedure (usually 1 or 2 hours).
4. Rinse with buffer 1 (3 X 10 mins).
5. Incubate with a 1/100 to 1/200 dilution of the UNDECAGOLD reagent in buffer 2 for 2 hours at room temperature.
6. Rinse with buffer 3 (3 X 5 mins), then buffer 4 (2 X 5 mins).
7. OPTIONAL (may improve sensitivity): Postfix with glutaraldehyde, 1 % in buffer 4 (10 mins).
8. Rinse with deionized water (2 X 5 mins).
9. OPTIONAL (may reduce background): Rinse with 0.05 M EDTA at pH 4.5 (5 mins).
10. Develop with freshly mixed silver developer for 20-25 minutes or as directed in the instructions for the silver reagent, twice. Rinse thoroughly with deionized water between developments to remove all the reagent.
11. Rinse several times with deionized water.

CAUTION: UNDECAGOLD particles degrade upon exposure to concentrated thiols such as β-mercaptoethanol or dithiothreitol. If such reagents must be used, concentrations should be kept below 1 mM and exposure restricted to 10 minutes or less.

Buffer 1: 20 mM phosphate
150 mM NaCl
pH 7.4
4% BSA (bovine serum albumin)
2 mM sodium azide (NaN₃)

1% normal serum; use serum of the host animal
for the UNDECAGOLD-conjugated
antibody

0.1% gelatin (Type B, approx. 60 bloom)

Optional, may reduce background:

0.5 M NaCl
0.05% Tween 20

Buffer 2: 20 mM phosphate
150 mM NaCl
pH 7.4
0.8% BSA

Buffer 3: 20 mM phosphate
150 mM NaCl
pH 7.4
0.8% BSA (bovine serum albumin)
2 mM sodium azide (NaN₃)

Buffer 4 (PBS):
20 mM phosphate
150 mM NaCl
pH 7.4

Other procedures may be used; for example the UNDECAGOLD reagent may be used as a tertiary labeled antibody, or a custom UNDECAGOLD conjugate may be the primary antibody. If additional antibody incubation steps are used, rinse with buffer 3 (3 X 10 mins) after incubation.

REFERENCES

1. Furuya, F. R., and Hainfeld, J. F.; *J. Histochem. Cytochem.*, **40**, 177 (1992); Furuya, F. R., Hainfeld, J. F., and Powell, R. D., *Proc. 49th Ann. Mtg., Electron Micros. Soc. Amer.*; G. W. Bailey and Hall, E. L. (Eds.); San Francisco Press, San Francisco, CA, p. 286 (1991).
2. Lipka, J. J., Hainfeld, J. F., and Wall, J. S., *J. Ultrastruct. Res.*, **84**, 120 (1983).
3. D'Souza, S. E., et al, *J. Biol. Chem.*, **263**, 3943 (1988).
4. J. E. Beesley, in "Colloidal Gold: Principles, Methods and Applications," M. A. Hayat, ed., Academic Press, New York, 1989; Vol. 1, pp 421-425.
5. Wall, J. S., Hainfeld, J. F., Bartlett, P. A., and Singer, S. J.: *Ultramicroscopy*, **8**, 397 (1982).
6. Moeremans, M. et al., *J. Immunol. Meth.* **74**, 353 (1984).

Technical Assistance Available.

For a complete list of references citing this product, please visit our world-wide-web site at <http://www.nanoprobes.com/Ref.html>.